

How Many Quantibody Slides Do I Need?

1. Running the full set of standards on every slide (Ideal)

of samples / 8

2. Running the full set of standards on every other slide (Recommended)

of samples / 11





3. Running the full set of standards on every 3 slides

of samples / 12

4. Running the full set of standards on every 4 slides (Maximum)

of samples / 12.5


Full Set of Standards Run on Every Slide (Ideal)

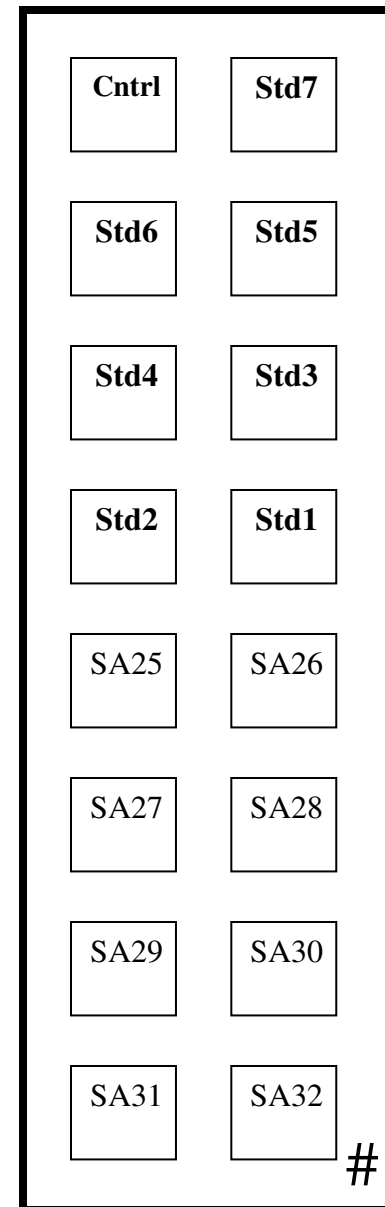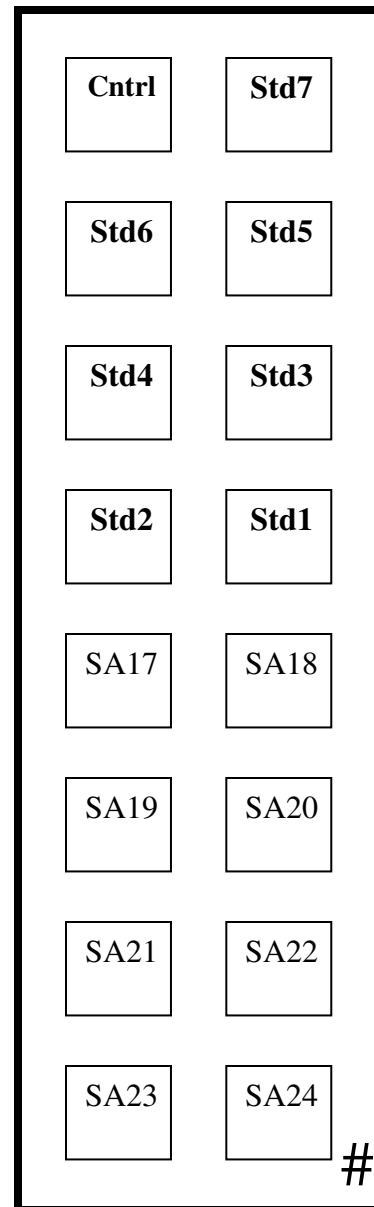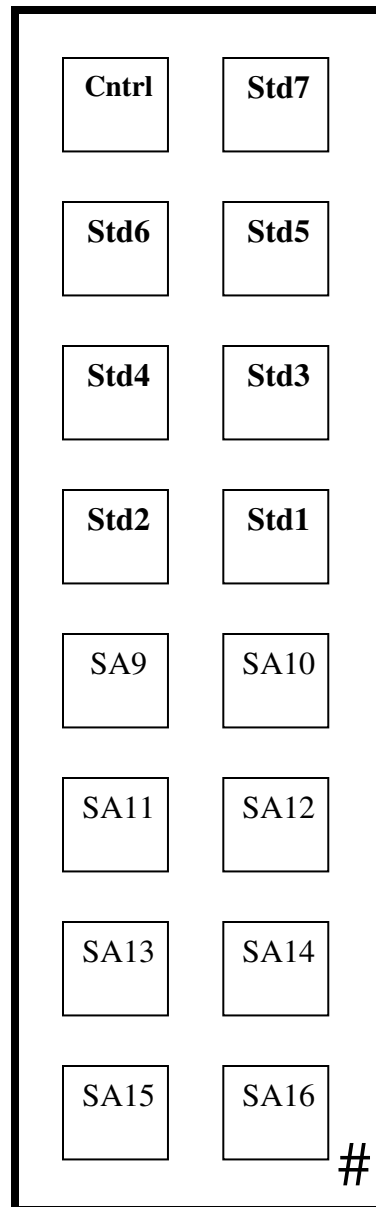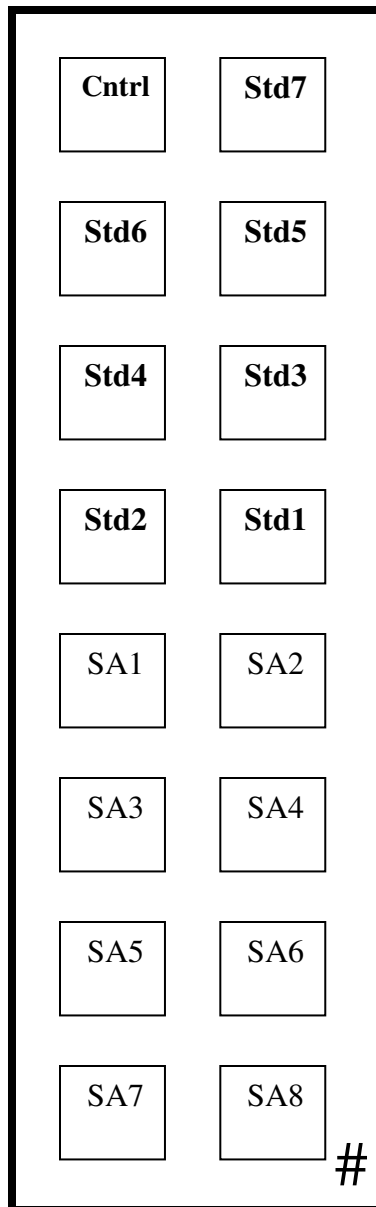

Full Set of Standards Run on Every 2 Slides (Recommended)

| | | | | | | | |
|-------|------|-------|------|-------|------|-------|------|
| Cntrl | Std7 | Cntrl | Std3 | Cntrl | Std7 | Cntrl | Std3 |
| Std6 | Std5 | SA9 | SA10 | Std6 | Std5 | SA31 | SA32 |
| Std4 | Std3 | SA11 | SA12 | Std4 | Std3 | SA33 | SA34 |
| Std2 | Std1 | SA13 | SA14 | Std2 | Std1 | SA35 | SA36 |
| SA1 | SA2 | SA15 | SA16 | SA23 | SA24 | SA37 | SA38 |
| SA3 | SA4 | SA17 | SA18 | SA25 | SA26 | SA39 | SA40 |
| SA5 | SA6 | SA19 | SA20 | SA27 | SA28 | SA41 | SA42 |
| SA7 | SA8 | SA21 | SA22 | SA29 | SA30 | SA43 | SA44 |
| | # | | # | | # | | # |

Full Set of Standards Run on Every 3 Slides

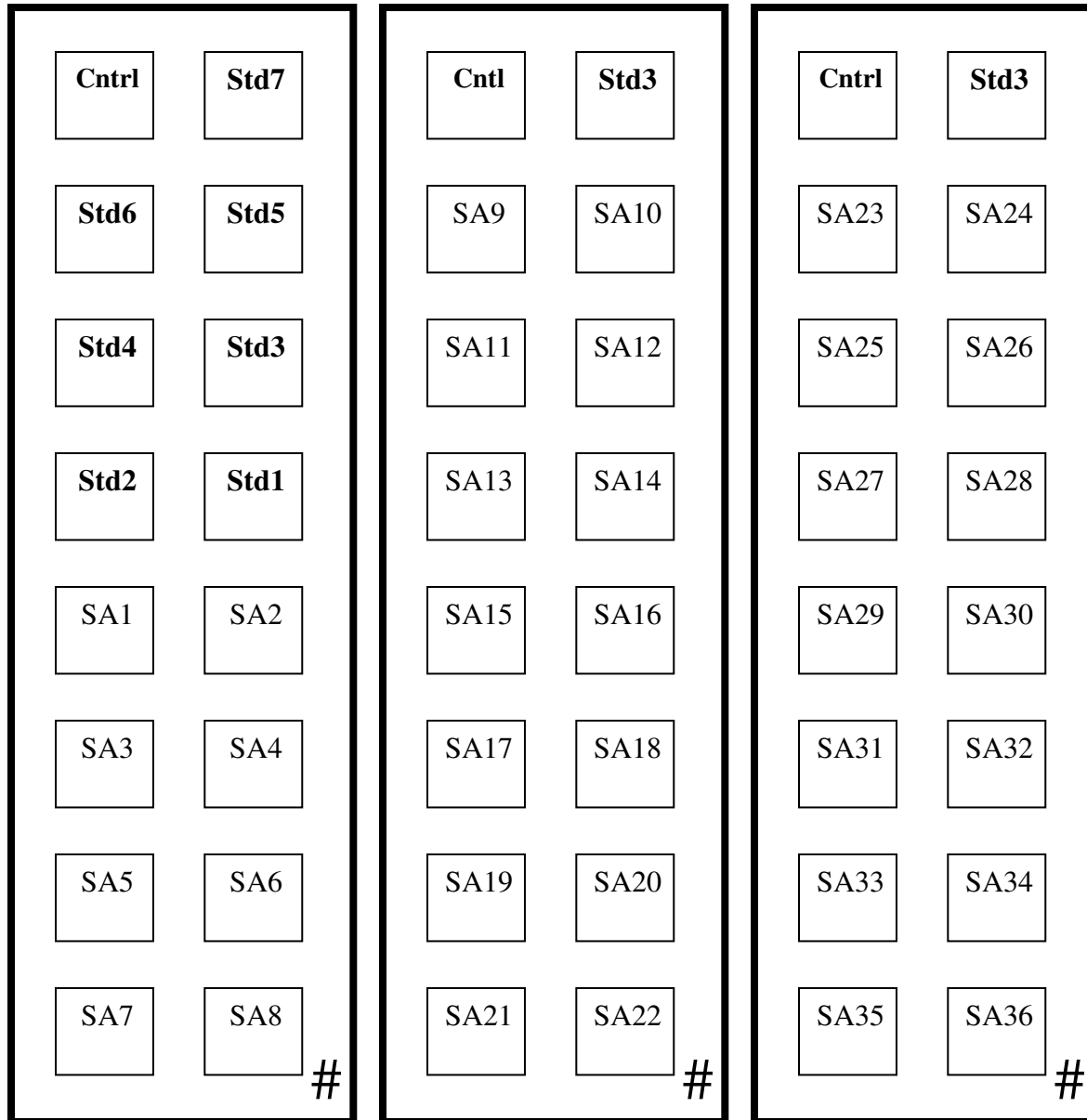

Full Set of Standards Run on Every 4 Slides (Maximum)

| | | | | | | | |
|-------|------|-------|------|-------|------|-------|------|
| Cntrl | Std7 | Cntrl | Std3 | Cntrl | Std3 | Cntrl | Std3 |
| Std6 | Std5 | SA9 | SA10 | SA23 | SA24 | SA37 | SA38 |
| Std4 | Std3 | SA11 | SA12 | SA25 | SA26 | SA39 | SA40 |
| Std2 | Std1 | SA13 | SA14 | SA27 | SA28 | SA41 | SA42 |
| SA1 | SA2 | SA15 | SA16 | SA29 | SA30 | SA43 | SA44 |
| SA3 | SA4 | SA17 | SA18 | SA31 | SA32 | SA45 | SA46 |
| SA5 | SA6 | SA19 | SA20 | SA33 | SA34 | SA47 | SA48 |
| SA7 | SA8 | SA21 | SA22 | SA35 | SA36 | SA49 | SA50 |
| | # | | # | | # | | # |